

## **ANTIGENIX AMERICA Inc.**

### Product Information Summary

#### **EZPrep Cell Strainer (35 micron)**

**DESCRIPTION:** Prepare Cell Suspensions for Flow Cytometry Analysis.

Cat No: EZP35S            Qty: 50    \$95.00

Cat No. EZP35L           Qty: 250   \$350.00

**APPLICATION:** The EZPrep Cell Strainer, 35 micron, can be used to remove cell aggregates or particles that are larger than 35  $\mu\text{m}$  in cell suspension preparations for flow cytometry analysis.

#### **PROCEDURE:**

1. Place an EZPrep Cell Strainer in a 12x75mm test tube.
2. If needed, seat the filter disc using the open end of a P1000 pipette tip.
3. Add up to 2 mL of cell suspension.
4. Cover the open end with gloved thumb to seal the strainer.
5. Gently squeeze the strainer to push the sample into the test tube
6. The flow-through fraction, with particles smaller than 35  $\mu\text{m}$ , is ready for flow cytometry analysis



**STORE:** Room Temp.    Research Use Only    [WWW.ANTIGENIX.com](http://WWW.ANTIGENIX.com)